



Transfection Efficiency of NeuroFECT™ Reagent on Rat Hippocampal Cells

A. NeuroFECT reagent only

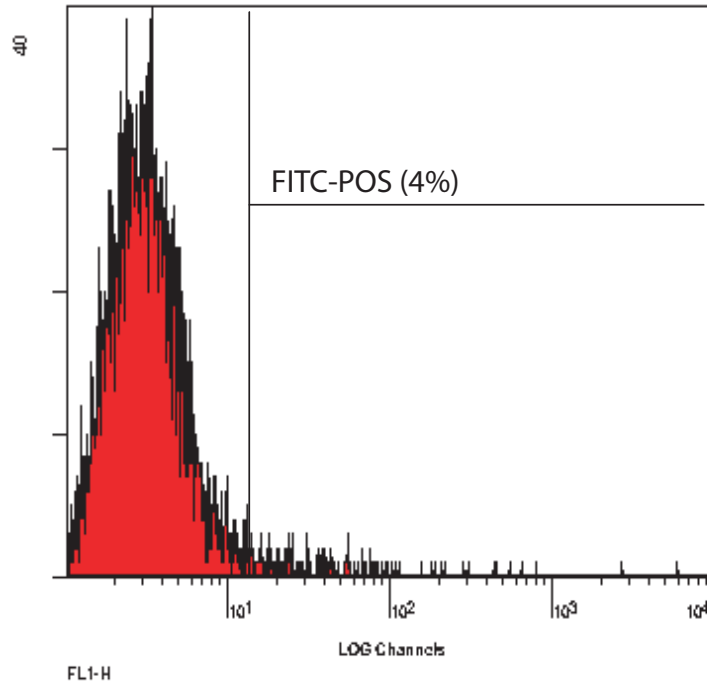

B. NeuroFECT reagent and gWiz-GFP expression vector

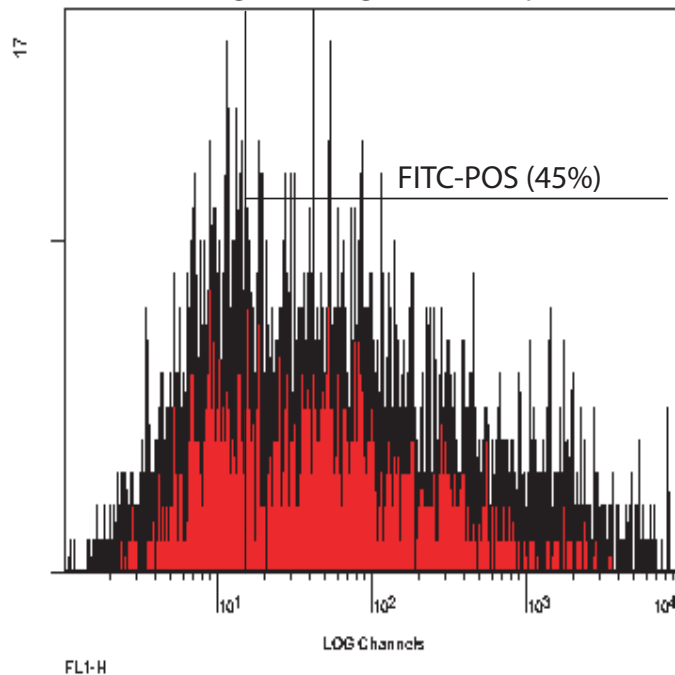

Primary E18 rat hippocampal cells were transfected with 0.7 μ l NeuroFECT and 0.5 μ g gWiz-GFP plasmid. Transfection efficiency was measured by FACS analysis 48 hours post-transfection. Graph B shows a 45% transfection efficiency.



10190 Telesis Court, San Diego, CA 92126
Toll-Free 888-428-0558 or 858-457-1919
Fax 858-623-9494 or 858-558-3617
genlantis.com