



# Same-Day Transfection using GenePORTER® 3000 Transfection Reagent


1. Plate log-phase cells in complete medium and incubate at 37°C in 5% CO<sub>2</sub> for 30 minutes.

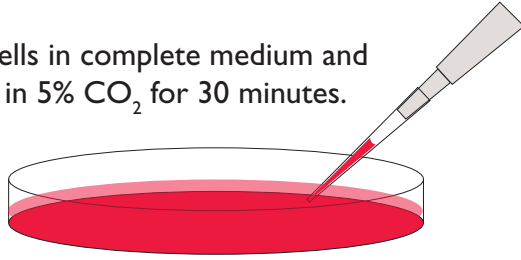


2. During incubation, prepare GenePORTER 3000 lipoplexes.

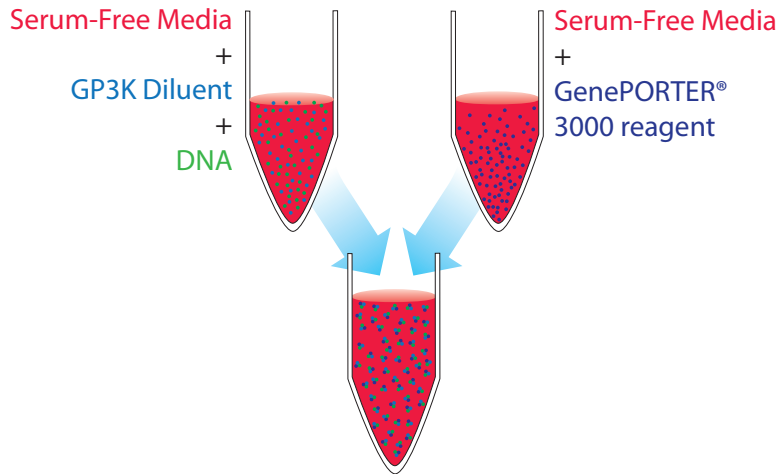


3. Remove complete medium. Replace with a equal volume of serum-free medium. Add GP3K lipoplexes.

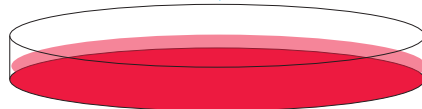

4. Incubate cells for 4 hours

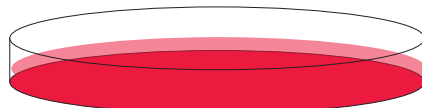

5. Add equal transfection volume complete medium with 20% serum.

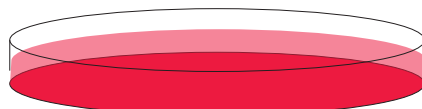

6. Assay after 24-72 hours.

