

cf-DNA/cf-RNA Preservation Tube Appearance at Various Stages

Before collection


After collection

After 5X Inversion

Resting 1 hour

After RT Shipping

After Centrifugation

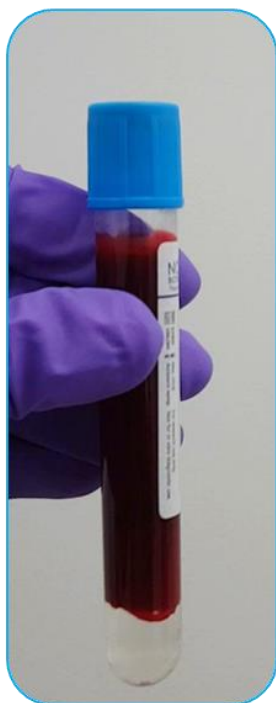

Note: Unlike other collection tubes blood will immediately start to settle and condense in the bottom of the tube leaving a cell-free supernatant. This is completely normal with the Norgen chemistry and does not indicate clotting. Following 5X inversion to mix and left to settle the dense lower fraction will not mix. Simply centrifuge as in the protocol and proceed with isolation of cf-DNA or cf-RNA from plasma.